

## Santai Adaptor Kit for 800g , 1600g and 3kg Flash Columns

### Procedure –

1. Luer connector kit for large columns (800g, 1600g and 3kg) (Santai Item#SLL-0102 and BLL-0102) consists of four main components as follows.



← Luer Connector for Inlet



← Luer Connector for Outlet

¼-28 Female to Male Luer →





¼-28 Female to Female Luer →



Namely

- Luer Connector for Outlet
- Luer Connector for Inlet
- ¼-28 Female to Male Luer
- ¼-28 Female to Female Luer

Product Picture	Santai Item#	Description
	SLL-0102	Luer connector kit for standard columns (4g-330g)
	BLL-0102	Luer connector kit for large columns (800g, 1600g and 3kg)

2. Install the Luer connector for outlet at the outlet of the column. Install the ¼-28 Female to Female Luer on it. Please refer the below pictures.



3. Install the Luer connector for inlet at the inlet of the column. This can be installed by hand and needs to be screwed in manually. Install the ¼-28 Female to male Luer.



4. The entire assembly will look like as shown in the below picture. Connect the inlet and outlet tubings from the Flash system to the outlet and inlet of the column.

